

Cellular-resolution Photostimulation



Polygon

DMD Pattern Illuminators

INTRODUCTION

The Polygon DMD pattern illuminators are Mightex's market-leading modules for targeted photostimulation. The Polygon provides precise spatio-temporal control of light with subcellular resolution, making it the perfect illumination tool for scientific research. Compatible with any upright or inverted microscope, the Polygon enables researchers to send light to anywhere on their specimen, and in any shape, size and complexity now within a large projection field of view. In addition, multiple regions-of-interest (ROIs) can be illuminated simultaneously, and patterns can be switched at kHz speeds. Different wavelengths of light can be used with the Polygon for virtual simultaneous multi-color illumination of unique ROIs. Polygon systems seamlessly integrate via TTL with other equipment such as electrophysiology tools or cameras.

DMD TECHNOLOGY

The Polygon uses digital micromirror device (DMD) technology to illuminate multiple ROIs simultaneously. A DMD chip is composed of up

to millions of micromirrors that can be individually turned ON/OFF to reflect light onto the sample. Thus, you can assign each mirror to control the area(s) of illumination and create any number of different-sized ROIs simultaneously.

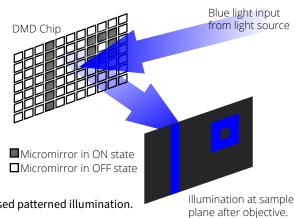


Figure 1. DMD-based patterned illumination.

FEATURES

- Illuminate any Shape or Size Within Large Field of View
- O Multi-Wavelength Illumination of **Distinct ROIs**
- O Simultaneous Multi-Region Illumination
- Fast Pattern Switching Speed
- Fits on any Microscope
- External Equipment Synchronization

APPLICATIONS

- O Neuroscience: Single-cell **Resolution Optogenetics**
- O Cell Biology: Subcellular **Resolution Optogenetics**
- Freely-Behaving Optogenetics
- Cortex-Wide Optogenetics
- Photoactivation. Photoconversion & Photoswitching
- Uncaging
- Photopatterning






POLYGON1000

MORE PIXELS, FASTER SPEED SAME VERSATILITY

Larger field of view. Finer resolution.

> New, larger DMD chip combined with front tube optics enables larger field of view without compromising resolution and power.

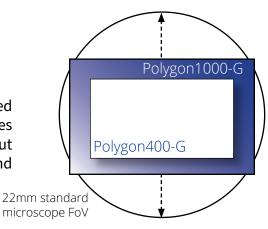

82% increase in illumination field of view*

Finer pixel resolution within larger field of view*

*Comparison between Polygon400-G model and large field-of-view Polygon1000-G.

Faster than anything else in the market.

Increased maximum frame rate means better temporal resolution for advanced physiologically-relevant experiments and virtually simultaneous 2-color illumination of distinct ROIs.

1000 SERIES

6,600 fps

400 **S**ERIES

4,000 fps

1.65X faster pattern switching speeds

Real-time projection. Closed-loop experiments.


> Faster uploading time enables the Polygon1000 to perform real-time pattern illumination for closed-loop experiments (faster than HDMI refresh rates [60Hz/120Hz]).





More power. Extra flexibility.

Larger chip and improved optics enhance transmission efficiency enabling the Polygon1000 to achieve increased power density at the specimen level, and giving the researcher more room for intensity control.

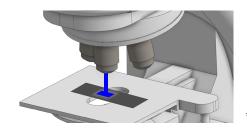



*Comparison between Polygon400-G model and Polygon1000-G.





POLYGON MODELS

1000 SERIES

POLYGON1000-G

P/N: DSI-K3-000

- Accepts a 3mm-core liquid lightguide.
- Can be used with any light source.

- Wavelength range: 350-700nm.
- Add-on front tube available for large field of view.

POLYGON1000-DL

P/N: DSI-K3-L20

- Accepts SMA-connectorized optical fiber patch cord Wavelength range: 400-700nm. (400µm, 0.22NA recommended).
- Compatible with laser sources.





400 SERIES§

POLYGON400-G

P/N: DSI-G-YYY



- Accepts a 3mm-core liquid lightguide.
- Can be used with any light source.
- Wavelength range: 400-700nm.






MICROSCOPE INTEGRATION

The Polygon can be coupled to most commercially available inverted and upright microscopes (Nikon, Leica, Zeiss, Olympus) in the following configurations:


INFINITY PATH CONFIGURATION

This configuration projects the spatial patterns at infinity, and hence it is mounted directly into the infinity path of a microscope by using a beam-combiner (for upright microscopes) along with an adaptor that matches the exact make/model of the microscope.

UPRIGHT

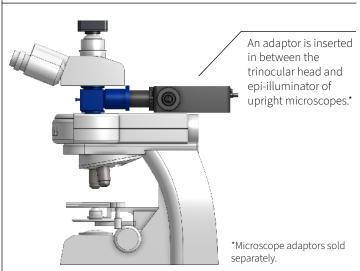

INVERTED

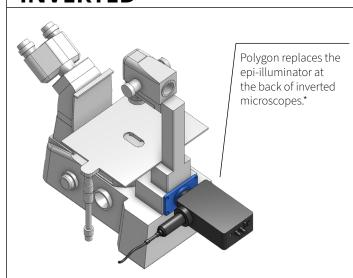

PLEASE CONTACT MIGHTEX FOR INTEGRATION OF MULTIPLE POLYGONS AND ALTERNATIVE MICROSCOPE INTEGRATION SOLUTIONS.

IPX INFINITY PORT EXPANDER

Mightex's IPX expands an infinity-path port on a microscope into a maximum of 4 ports. It is compatible with all Mightex Polygon models as well as with Mightex and 3rd party widefield epi-fluorescent illumination sources via standard 3mm core liquid lightguide. It also supports cameras and laser scanners via appropriate adaptors. Ports 2, 3 and 4 feature pitch-yaw adjustable dichroic holders for centering FOV and each port can be mounted on either sides of the main IPX chassis, to avoid mechanical conflict with surrounding environment.

C-MOUNT CONFIGURATION

If the infinity path of your microscope is unavailable, this configuration can be mounted onto one of the standard C-mount camera ports of your microscope.



LAPP CONFIGURATION

Do you have a Nikon microscope with a LAPP modular illumination system? We provide a Polygon format that is LAPP compatible. Please contact Mightex for more information.





TECHNICAL SPECIFICATIONS

| ILLUMINATION FIELD-OF-VIEW & RESOLUTION | | | | | | | | |
|---|--------------------|-------------------------------|---|-------|---------|-------|--|--|
| Model | Field of View | Projection Area Dimensions | Commercial Microscope (1X Objective) ^a | | | | | |
| | | | Leica | Nikon | Olympus | Zeiss | | |
| 1000 SERIES | | | | | | | | |
| POLYGON1000-G | Standard | Height mm | 6.2 | 6.2 | 5.5 | 5.1 | | |
| | | Width mm | 9.9 | 9.9 | 8.9 | 8.1 | | |
| | | Diagonal mm | 11.6 | 11.6 | 10.5 | 9.6 | | |
| | | Pixel Size μm | 7.6 | 7.6 | 6.9 | 6.3 | | |
| | Large ^c | Height mm | 12.4 | 12.4 | 11 | 10.2 | | |
| | | Width mm | 19.8 | 19.8 | 17.8 | 16.2 | | |
| | | Diagonal mm | 23.2 | 23.2 | 21 | 19.2 | | |
| | | Pixel Size μm | 15.2 | 15.2 | 13.8 | 12.6 | | |
| POLYGON1000-DLb | Standard | Diameter ^b mm | 12.4 | 12.4 | 11 | 10.2 | | |
| | | Pixel Size μm | 15.2 | 15.2 | 13.8 | 12.6 | | |
| 400 SERIES [§] | | | | | | | | |
| POLYGON400-G | Standard | Height mm | 8.7 | 8.7 | 7.8 | 7.2 | | |
| | | Width mm | 15.5 | 15.5 | 13.9 | 12.7 | | |
| | | Diagonal mm | 17.7 | 17.7 | 16.0 | 14.6 | | |
| | | Pixel Size μm | 18.0 | 18.0 | 16.2 | 14.8 | | |

^a To calculate illumination field-of-view and pixel resolution at the specimen, simply divide the above numbers by the magnification of the objective.

| CONTROL & TIMING | | | | | | |
|---------------------------------------|---------------------------------|-------------|--|--|--|--|
| | 400 SERIES | 1000 SERIES | | | | |
| Maximum Frame Rate ^a fps | 4,000 | 6,600 | | | | |
| Input Trigger | TTL, BNC connector | | | | | |
| Input Trigger Current | Minimum 8mA | | | | | |
| Input Trigger Delay μs | 50 | | | | | |
| Output Trigger | tput Trigger TTL, BNC connector | | | | | |
| Output Trigger Delay | User Programmable | | | | | |
| Input Uploading Speed ms/frame | 150 | 4 | | | | |

 $^{^{\}rm a}$ Values at 1bit depth. For grayscale features please contact Mightex for more information.



^b Polygon1000-DL has a circular illumination field-of-view.

^c Large field-of-view front tube lens required. Sold separately.

[§] Polygon400 series has been discontinued and replaced with the 1000 Series.



SYSTEM & COMMUNICATION

| Operating System§ | Windows XP, Vista, 7, 8, and 10 | | | | |
|-------------------|---|--|--|--|--|
| Software | Mightex PolyScan2 Nikon NIS Elements NeuroPG* | | | | |
| | SuperSpeed USB3.0 | | | | |
| Power Supply | 5Vdc 3A input power | | | | |
| Screen Resolution | 1,366x768 or higher | | | | |

[§] Polygon1000 supported by 64bit systems only.

ORDER NOW

Our primary goal is to help you find the optimal solution for your research. We have a dedicated technical support and sales team committed to providing expert guidance on our Polygon models and other Mightex products.



Please visit www.mightexbio.com/contact to submit an inquiry form today!

CONTACT US

US OFFICE

1241 Quarry Lane, Suite 105 Pleasanton, CA 94566 **USA**

TEL: 1-925-218-1885

CANADA OFFICE

111 Railside Road, Suite 201 Toronto, ON M3A 1B2 Canada

TEL: 1-416-840-4991





www.mightexbio.com

^{*}Polygon400-G only